

Foots installation



一、Foots installation

Required Tools:

- 1.No.36 Wrench (picture 1)
- 2.Foots: 4 (picture 2)





二、Operating steps:

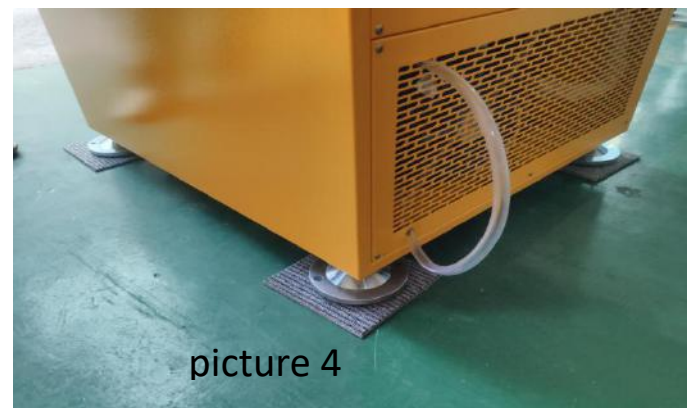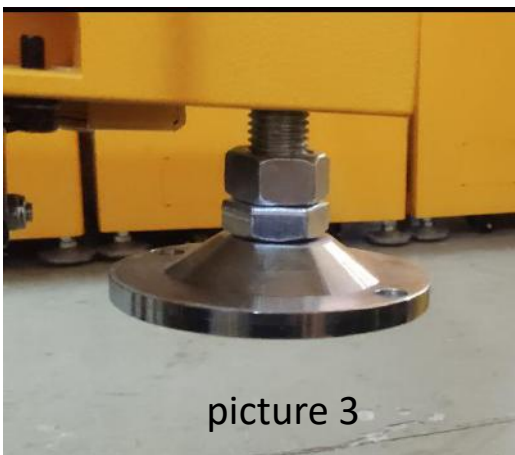

1. There are two screws, the lower one is used to adjust the height, the upper one is used to fix. (picture 3)
2. Put the foot pads under the four feet. (picture 4)

3. The lower screw is turned in below clockwise by 36 wrench until the caster can leave the ground .When tightening the foot, hold the wrench in one hand and the caster in the other hand. (picture 5,picture 6)

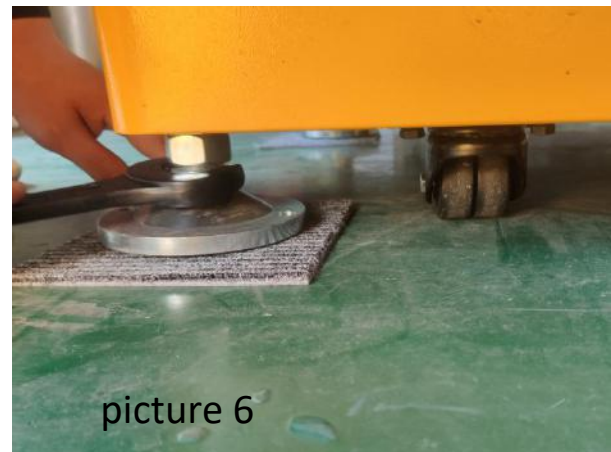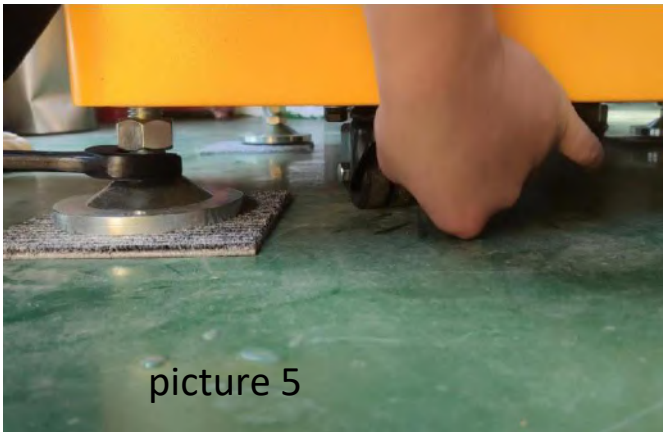

4. If it has 3 units, first to turn on the middle and bottom units and adjust the speed to 180rpm. Then you can touch the feet by your hand and make small adjustments when you feel shaking.



5. when you don't feel shaking about four feet, turn the upper screw. (Picture 8)
(Note: Do not tighten at this time, reserve some space that can be loosened.)
6. Turn on the third unit and adjust the speed to 180 rpm. Then you can touch the feet by your hand and make small adjustments when you feel shaking.
7. When the 3 units are running at the same time and the foots are not shaking, the upper screw can be tightened. (Picture 9)

