

粘板的清洗方法与注意事项

Sticky mat cleaning methods and precautions

1. 防止紫外线照射（紫外照射降低使用寿命），如需使用紫外灯，必须贴好粘板原保护膜。

1. Protect from UV light (UV exposure will effect service life). If you need to use the UV lamp, the original protective film of the sticky mat must be pasted.

2. 清洗粘板：建议使用超声波清洗（如加洗洁精，请选用中性洗洁精清洗），自然晾干即可。

2.Cleaning the sticky mat: It is recommended to clean with ultrasonic (if detergent is added, please use neutral detergent for cleaning), then dry naturally.

3. 禁止使用布或纸巾等物品擦拭。

3.Do not use cloth or paper towels to wipe.

4. 切勿用酒精等用品消毒或清洗。

4. Do not disinfect or clean with items such as alcohol.